

SNV/U High Performances UV Microchip Series



KEY FEATURES

- 355 nm and 266 nm
- Repetition rate up to 20 kHz
- Ultrashort pulses down to 550 ps
- Multi-kW peak power
- Excellent beam quality
- Efficient, air-cooled
- Sealed package, extremely long life

For generating high peak power ultraviolet pulses of a few hundred picoseconds, microchip lasers are economical, compact, and reliable. Micro-joule UV pulses are generated by harmonic conversion of the IR passively Q-switched Nd:YAG engine. Microchips are also easy to operate and service ; controllers can be used with every laser head model and swapped within minutes while conserving constant performances. The SNV and SNU series are designed for high average power, delivering multi-kW peak power at repetition rates up to 20 kHz.

APPLICATIONS

- Semiconductor inspection
- Laser-induced fluorescence (LIF)
- Micro dissection
- Organic compound marking and micromachining

- Biohazard detection
- Time resolved fluorescence
- Laser Induced Breakdown Spectroscopy (LIBS)



TECHNICAL SPECIFICATIONS

	SNV-05P-100	SNV-20F-100 ⁽⁷⁾	SNU-02P-100	SNU-20F-100
Wavelength	355nm	355nm	266nm	266nm
Repetition Rate	>5kHz	>19kHz	>6kHz	>19kHz
Constant Pulse width range (FWHM) ⁽¹⁾	<0.6ns	<0.6ns	<0.6ns	<0.6ns
Output power ⁽²⁾	>5mW	>10mW	>2mW	>10mW
Output energy	>0.5µJ	>0.5µJ	>0.3µJ	>0.5µJ
Peak Power	>0.7kW	0.7kW	>0.5kW	>0.7kW
Short term (1min) power stability ⁽³⁾	<±1%	<±1%	<±1%	< <u>±</u> 2%
Long term (6 hrs) power stability ⁽³⁾	<±5%	<±5%	<±5%	<±5%
Beam profile	Gaussian TEM00	Gaussian TEM00	See note (5)	See note (5)
Full angle divergence Horizontal@1/e ² Vertical@1/e ²	8.5±2mrad 6±2mrad	11±2mrad 7±2mrad	11±2mrad <1.5mm ⁽⁶⁾	11.5±2mrad 0.65±0.25mrad
M ²⁽⁴⁾	<1.3	<1.3	<1.3	<1.4
Gaussian fit in far field	N/A	N/A	N/A	>85%
Polarization	Linear PER>20dB	Linear PER>20dB	Linear PER>20dB	Linear PER>20dB
Package dimensions	180x55x36mm	186x60x36mm	180x55x36mm	210x60x36mm
Package weight	400g	500g	400g	500g
Options (table p3)	С	С	С	С
Options included	-	S	-	S

NOTES

(1) Measured with 1Ghz photodiode and 1GHz/10GS/s oscilloscope.

(1) measured with 1Ghz photodiode and 1GHz/10GS/s oscilloscope. (2) Measurement performed with an OPHIR thermal power sensor (OPHIR 3A-FS-SH) (3) For temperature variation < \pm 3°C and < 3°C/hour, stability is measured with calorimeter - detector band [DC, 2Hz] (4) Mean average value M = $\sqrt{(XY)}$, X and Y being respectively the major and minor axis of the ellipse (5) Beam exhibits different profile in horizontal (Gaussian) and vertical ((sin x /x)² in far-field) plan (6) 5%/95% diameter, at 300mm from laser output (7) Contact factory for availability



COMPLEMENTARY INFORMATION & OPTIONS

Environment Parameters				
Operating Temperature Range	15-35°C			
Maximum Laser Head Baseplate Temperature	<50°C			
Maximum Power Consumption	<40W			
Laser Head Thermal Dissipation	<15W			
Storage Temperature	0-50°C			
Shock of 11ms according to IEC 68-2-27, non operating	25g			
Vibration 5Hz to 500Hz sinusoïdal according to IEC 68-2-6	2g			

Certification					
Laser classification according to IEC 60825-1:2007	3B for SNV-05P and SNV-20F 4 for SNU-02P and SNU-20F				
CDRH	Yes, if used with a -DR1 controller				
ROHs	Yes				


Options					
Collimation (C)	With collimated beam				
Synchronization output (S)	TTL compatible output signal for synchronization/monitoring				

Available Controller Types					
Model	Туре	Input Power	CDRH		
MLC-03A-DR1	Desktop	100-240 V AC	Yes		
MLC-03A-MR1	Module	12 V DC	No		
MLC-03A-BR1	Board	12 V DC	No		

All information contained herein is believed to be accurate and is subject to change without notice. Teem Photonics may not be held liable for its use. Teem Photonics, its subsidiaries and affiliates, reserve the right to modify or withdraw, at any time and without any notice, specifications, product design, product component. Some options may not be available for all products. Please contact Teem Photonics for details. August 18



CDRH LASER HEAD MECHANICAL DRAWINGS: SNV-05P-100

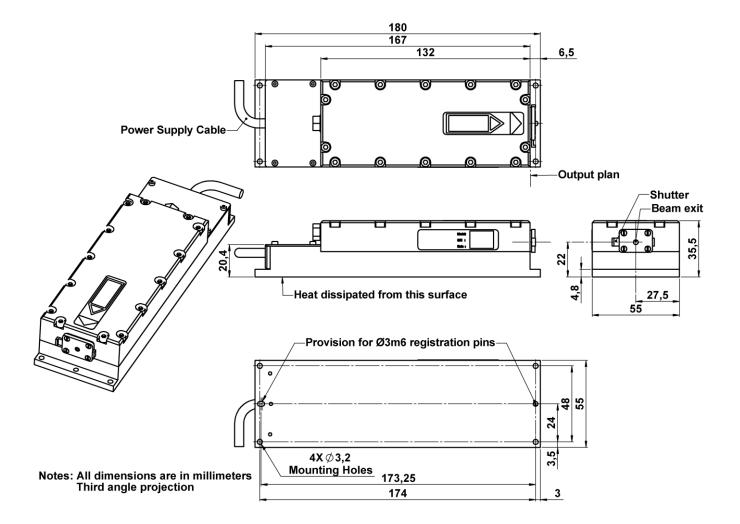


All information contained herein is believed to be accurate and is subject to change without notice. Teem Photonics may not be held liable for its use. Teem Photonics, its subsidiaries and affiliates, reserve the right to modify or withdraw, at any time and without any notice, specifications, product design, product component. Some options may not be available for all products. Please contact Teem Photonics for details. August 18

www.teemphotonics.com



CDRH LASER HEAD MECHANICAL DRAWINGS: SNU-02P-100

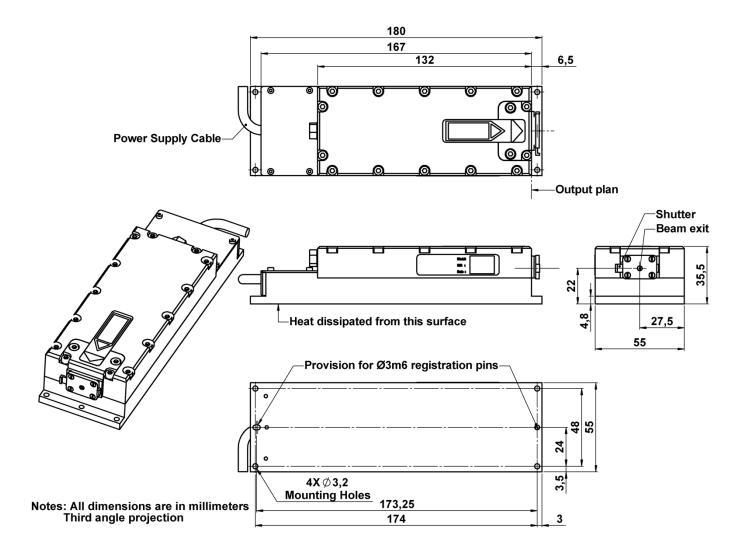


All information contained herein is believed to be accurate and is subject to change without notice. Teem Photonics may not be held liable for its use. Teem Photonics, its subsidiaries and affiliates, reserve the right to modify or withdraw, at any time and without any notice, specifications, product design, product component. Some options may not be available for all products. Please contact Teem Photonics for details. August 18

www.teemphotonics.com SNV & SNU I 5/7



CDRH LASER HEAD MECHANICAL DRAWINGS: SNV-20F-100

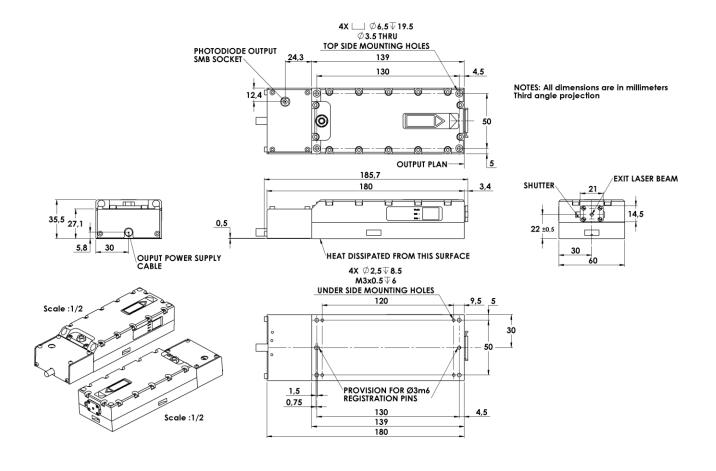


All information contained herein is believed to be accurate and is subject to change without notice. Teem Photonics nay not be held liable for its use. Teem Photonics, its subsidiaries and affiliates, reserve the right to modify or vithdraw, at any time and without any notice, specifications, product design, product component. Some options may to be available for all products. Please contact Teem Photonics for details. August 18

www.teemphotonics.com SNV & SNU I 6/7



CDRH LASER HEAD MECHANICAL DRAWINGS: SNU-20F-100

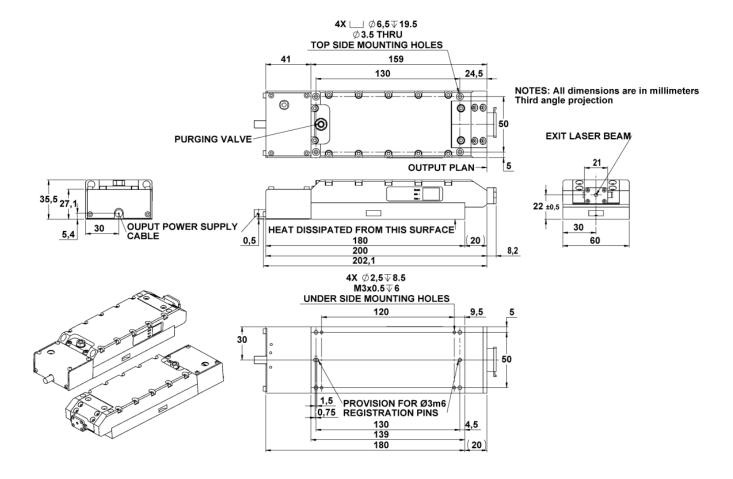

All information contained herein is believed to be accurate and is subject to change without notice. Teem Photonics may not be held liable for its use. Teem Photonics, its subsidiaries and affiliates, reserve the right to modify or withdraw, at any time and without any notice, specifications, product design, product component. Some options may not be available for all products. Please contact Teem Photonics for details. August 18

www.teemphotonics.com SNV & SNU I 7/7